

## Ettan DIGE Imager

The Ettan™ DIGE Imager (Fig 1) is a scanning CCD camera designed for applications in the life sciences. In particular, it has been engineered to create high quality images of 2-D DIGE gels. By combining high resolution with precise motion control, the Ettan DIGE Imager produces accurate multichannel images of your Cy™2-, Cy3-, and Cy5-labeled gel. The system has also been designed to image a wide range of other fluorescent gel applications, and can be used with the new ECL Plex™ Western Blotting System.

The imager is controlled from the Ettan DIGE Imager software, and can be set up for a variety of gel formats. Results from the Ettan DIGE Imager are directly compatible with ImageQuant™ TL, ImageMaster™ 2D Platinum, and DeCyder™ 2-D Differential Analysis (DeCyder 2-D) Software.

ImageMaster 2D Platinum v6.0 comes in two flavors: DIGE and non-DIGE. The newly developed DIGE-enabled version of the software comprises powerful and versatile tools for the analysis of images from DIGE experiments. Together with the Ettan DIGE Imager, ImageMaster 2D Platinum v6.0, DeCyder 2-D v6.5, and DeCyder Extended Data Analysis (EDA) Software create a powerful platform of image capture and analysis tools that offers flexibility, versatility, and adaptability for protein profiling.



Fig 1. Ettan DIGE Imager

### Ettan DIGE Imager features

- High sensitivity for imaging faint spots
- Software-selectable wavelengths
- Filter bandwidth optimized to minimize crosstalk
- Adjustable exposure level to optimize sensitivity
- Fast scanning for increased throughput
- Full 16 bits/pixel depth for accurate quantitation
- Sealed environment for scanning and protecting wet samples from drying
- Dedicated removable cassettes for handling and scanning gels—cassettes are available for naked gels or membranes and DALT and SE 600 series gel sandwiches
- Compatibility with Ettan Spot Picker
- Flat field calibration
- Compatible with image analysis software, such as ImageMaster, ImageQuant, or DeCyder software for quantitative evaluation of samples
- Two 1-GB/sec Ethernet connections for data transfer




## Principle of operation

Upon excitation by the metal halide lamp, light is emitted from a fluorescently labeled sample in proportion to the amount of labeled compound in the sample. Emitted light is collected and converted to an electronic signal in a CCD. The signal is displayed and analyzed. Data is stored in a 16-bit



TIFF to provide the digital resolution required to characterize subtle signal intensity differences over the wide dynamic range of the instrument.

Figure 2 shows a typical gel image produced with the Ettan DIGE Imager.

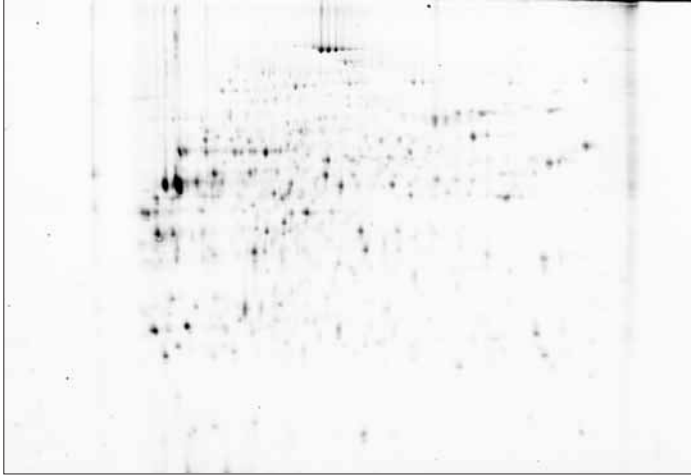

**Fig 2.** Cy2 image of a 2-D DIGE gel.

Figure 3 compares color images of the same 2-D DIGE gel acquired on the Ettan DIGE Imager and the Typhoon™ scanner. For 2-D DIGE gels, the two instruments produce images of similar quality.

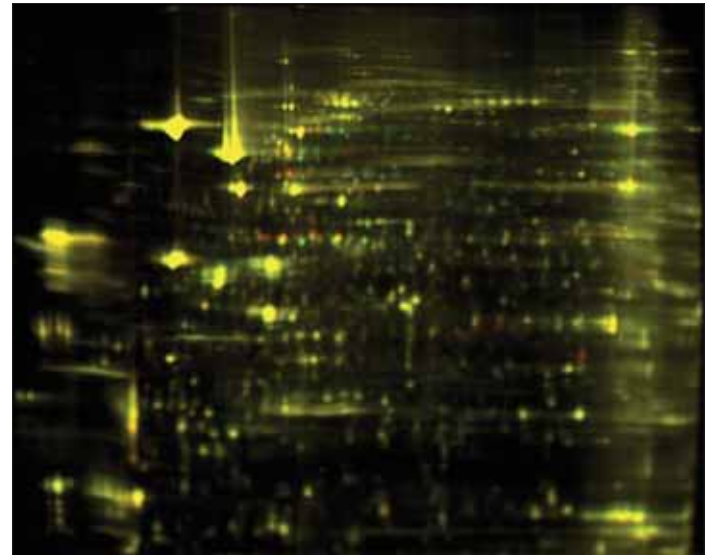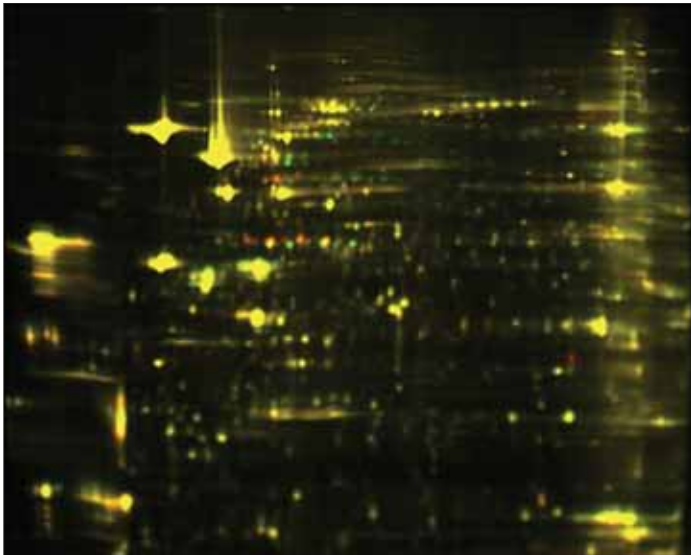

**Fig 3.** Images of a 2-D DIGE gel acquired on the Ettan DIGE Imager (left) and Typhoon scanner (right).

## Ettan DIGE Imager specifications

Light source	Metal halide lamp, coupled into fiber bundle
Lamp lifetime	1000 h
Scanning	Camera lens, moving sample, 1 × 1 cm images detected on-the-fly
Scanning time: 255 × 196 mm gel, three colors Ettan DALT gel, one color	Approximately 30 min 7–11 min
Scan area	270 × 205 mm
Uniformity	± 5% over entire scan area
Detection	Peltier-cooled CCD camera
Wavelength selection	Bandpass excitation (380–650 nm) and emission (465–695 nm) filters
Fluorescence dye compatibility	Cy2, Cy3, Cy5, Deep Purple™ Total Protein Stain and SYPRO™ Ruby
Detection limit	Cy2: 3 fmol; Cy3, Cy5: 1 fmol; Deep Purple Total Protein Stain: 1 ng; SYPRO Ruby: 5 ng
Linear detection range	3.5 orders of magnitude
Linearity	< 7.5% relative standard deviation over entire dynamic range
Pixel size	Software selectable: 40, 100, or 200 µm
Pixel accuracy	± 100 µm
Resolution	40 µm
Local reproducibility	± 50 µm between the first and third scans of the same gel
Signal reproducibility	< 5% difference between the first and third scans of the same gel
Software	Scan control software for Windows XP
External interface	10 Base-T Ethernet using TCP/IP protocol
Data format	16-bit (65,536 levels) TIFF, .GEL file extension
Instrument size	880 × 450 × 730 mm, footprint 710 × 680 mm
Weight	64 kg
Power requirements	115/230 V (autoswitching), 50–60 Hz, < 600 W

## Ordering information

Ettan DIGE Imager, including installation kit (power cable, CD with scan control software, test slides, and user manual with installation guidelines)	63-0056-42	<b>Imaging software</b>	
		ImageMaster 2D Platinum v6.0 DIGE Enabled	11-0034-25
		ImageMaster 2D Platinum v6.0	11-0034-27
Ettan DIGE Imager Cassette, 276 × 212 mm, for DALT gel sandwiches	11-0027-04	DeCyder 2-D Differential Analysis Software v6.5, preinstalled network (including PC and single concurrent network user license)	28-4012-01
Ettan DIGE Imager Cassette, 180 × 160 mm, for SE 600 series gel sandwiches	11-0027-32	DeCyder 2-D Differential Analysis Software v6.5, one network user license	11-0035-82
Ettan DIGE Imager Cassette, with low-fluorescent glass, for naked gels	11-0027-33	DeCyder Extended Data Analysis Software, one network user license	28-4012-03
		ImageQuant TL	63-0050-72

<b>Asia Pacific</b>	Tel: +852 2811 8693	Fax: +852 2811 5251
<b>Australasia</b>	Tel: + 61 2 9899 0999	Fax: +61 2 9899 7511
<b>Austria</b>	Tel: 01/57606-1619	Fax: 01/57606-1627
<b>Belgium</b>	Tel: 0800 73 888	Fax: 03 272 1637
<b>Canada</b>	Tel: 1 800 463 5800	Fax: 1 800 567 1008
<b>Central, East, &amp; South East Europe</b>	Tel: +43 1 982 3826	Fax: +43 1 985 8327
<b>Denmark</b>	Tel: 45 16 2400	Fax: 45 16 2424
<b>Finland &amp; Baltics</b>	Tel: +358-(0)9-512 39 40	Fax: +358 (0)9 512 39 439
<b>France</b>	Tel: 01 6935 6700	Fax: 01 6941 9677
<b>Germany</b>	Tel: 0761/4903-490	Fax: 0761/4903-405
<b>Italy</b>	Tel: 02 27322 1	Fax: 02 27302 212
<b>Japan</b>	Tel: +81 3 5331 9336	Fax: +81 3 5331 9370

<b>Latin America</b>	Tel: +55 11 3933 7300	Fax: + 55 11 3933 7306
<b>Middle East &amp; Africa</b>	Tel: +30 210 9600 687	Fax: +30 210 9600 693
<b>Netherlands</b>	Tel: 0165 580 410	Fax: 0165 580 401
<b>Norway</b>	Tel: 815 65 555	Fax: 815 65 666
<b>Portugal</b>	Tel: 21 417 7035	Fax: 21 417 3184
<b>Russia &amp; other C.I.S. &amp; N.I.S</b>	Tel: +7 (095) 232 0250 +7 (095) 956 1137	Fax: +7 (095) 230 6377
<b>South East Asia</b>	Tel: 60 3 8024 2080	Fax: 60 3 8024 2090
<b>Spain</b>	Tel: 93 594 49 50	Fax: 93 594 49 55
<b>Sweden</b>	Tel: 018 612 1900	Fax: 018 612 1910
<b>Switzerland</b>	Tel: 0848 8028 12	Fax: 0848 8028 13
<b>UK</b>	Tel: 0800 616928	Fax: 0800 616927
<b>USA</b>	Tel: +1 800 526 3593	Fax: +1 877 295 8102

[www.amershambiosciences.com/dige](http://www.amershambiosciences.com/dige)

**Amersham Biosciences Corp**

800 Centennial Avenue  
 PO Box 1327  
 Piscataway, NJ  
 08855-1327  
 USA

General Electric Company reserves the right, subject to any regulatory approval if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation. Contact your GE Representative for the most current information. © 2005 General Electric Company—All rights reserved. GE and GE Monogram are trademarks of General Electric Company. Amersham Biosciences, Cy, DeCyder, Deep Purple, ECL Plex, Ettan, ImageMaster, and ImageQuant are trademarks of GE Healthcare companies. SYPRO is a trademark of Invitrogen.



imagination at work